



NIS-Elements: Installing the Lumenera Driver

This document specifies the install procedures for the NIS-Elements Documentation (D) package and the Lumenera Driver for the Infinity 1-2C and the 1-3C.

Note: Operation of the Driver is only possible with the purchase of the Lumenera Driver material number 97152 (which is included in material # 97161 and 97162).

Note: Installing Lumenera Software is not required. Lumenera drivers are installed in Step #3.

Note number: 0013
Creation Date: 9/2010
Date Modified:
Software Version: 3.2
Package: D Package Only
Module: Lumenera Driver

1) Install NIS-Elements Documentation Package.

- We do not recommend installing the version on the shipped DVD, as a more recent version may be available on the download site specified in Step #2.
- Please note that the Lumenera driver will not be available in the NIS-Elements install 'Camera Section.'

2) The Lumenera driver is available for both camera modules at the NIS-Elements Download Site:

<http://www.nissoftware.net/NikonSaleApplication/>

Once at this site, enter User Name and Password:

User Name: Your Name (John Deer)

Password: the 8 digit code on the purple USB HASP key



After entering the site, scroll to the bottom of the page and click on 'NII Modules, Drivers and Macros:'



[NII Modules, Drivers and Macros](#)

After clicking on the above link, select and save the Lumenera zip folder that MATCHES the install build that you are installed in Step #1.

For example, if you installed Build 685, download the driver with the Build 685 in its file name:

[Lumenera-Install-Build-616.zip](#)

[Lumenera-Install-Build-641.zip](#)

[Lumenera-Install-Build-685.zip](#) ←

3) After opening and saving the zip file, double click on the 'Setup.exe' file.

- This will load the driver into the correct file locations automatically.

4) Open NIS-Elements Documentation Package by double clicking on icon.

5) Select the Lumenera Driver in the drop down menu at Start Up:

